



VitroView™ IHC Hemotoxylin Counterstain Kit (300 tests)

SKU No. VB-6004

Description: This ready-to-used (RTU) product is designed to be used after completion of immunohistochemical staining. The staining color will be light blue in nuclei.

Kit Contents

RTU Hematoxylin Solution 15ml ×2

Storage

Store at room temperature.

Application

Immunohistochemistry: Tested

Recommended Assay Procedure for Immunohistochemistry:

1. After DAB staining, incubate sections with 3 drops of RTU hematoxylin solution for 1-2 minutes
2. Rinse in tap water 2×2 min.
3. Dehydrate through 75% ethanol for 2 min, 95% ethanol for 2 min, and 100% ethanol for 2x3min. Clear in xylene for 2×5min.
4. Coverslip with mounting medium.

Note: This product is intended for research purposes only. This product is **not** intended to be used for therapeutic or diagnostic purposes in humans or animals. Avoid contact with eyes, skin and clothing. Do not ingest. Wear gloves.

Precautions

Handle with care in chemical hood. Avoid contact with eyes, skin and clothing. Do not ingest. Wear gloves.