



VivroView™ Permanent Aqueous Mounting Medium

SKU#: VB-8005

Description

VivroView™ Permanent Aqueous Mounting Medium is specially designed for the permanent preservation of tissue sections and cell smears immunohistochemically stained with chromogens that cannot be dehydrated using organic solvents.

Advantages

- With high refractive index
- No exposure to the organic fumes
- Permanent storage of slides and high resolution of tissue specimens

Application

Compatible with AEC, DAB, Fast Red, Oil Red, BCIP/NBT, BCIP/INT and fluorescent dyes and phycobiliproteins.

Package Size

15ml/ squeeze bottle

Storage

Product is stable for about 12 months at room temperature.

Usage Instruction

- Place slides flat on paper towel on countertop with specimen facing up.
- Blot excess water from the slide without letting tissue specimen dry prior to mounting.
- Apply 2-3 drops of Permanent Aqueous Mounting Medium to completely cover the specimen with a thin layer.
- Place a cover slip on the liquid and let sit at room temperature for 45 minutes to 1 hour to allow to dry.

Precautions:

- Avoid contact with skin and eyes.
- Harmful if swallowed.
- Follow all Federal, State, and local regulations regarding disposal.

Disclaimer:

This user manual serves as a general guideline. Users should adapt procedures based on specific experimental requirements and equipment specifications.