



VitroView™ Von Kossa Calcium Stain Kit

SKU: VB-3043

Introduction

The VitroView™ Von Kossa Calcium Stain Kit is a reliable and easy-to-use solution for detecting calcium deposits in tissue samples. Designed for use in paraffin-embedded sections, undecalcified bone, and calcium-rich tissues, this kit provides rapid, reproducible staining that highlights calcium salts in black or brown-black, with nuclei counterstained red and cytoplasm pink.

Ideal for histology, pathology, and biomedical research, the VitroView™ Von Kossa kit enables high-contrast visualization of calcium deposition, supporting studies of bone development, mineralization disorders, and tissue calcification. With optimized reagents and clear instructions, this kit delivers consistent, high-quality results for both routine and advanced research applications.

Key Benefits:

- Sensitive and specific detection of calcium salts
- Compatible with paraffin sections and undecalcified tissues
- Quick and reproducible staining
- Optimized for both UV light and standard laboratory lighting

Kit Contents

VB-3043-1	Silver Nitrate Solution	100 ml
VB-3043-2	Sodium Thiosulfate Solution	100ml
VB-3043-3	Nuclear Fast Red Solution	100ml

Storage

Silver Nitrate Solution: Stored at 2–8°C. The others: Stored at room temperature.

Procedure

1. Deparaffinize in xylene I for 6 minutes and II for 6 minutes.
2. Rehydrate
 - 1) Ethanol 100% (2 minutes)
 - 2) Ethanol 100% (2 minutes)
 - 3) Ethanol 95% (2 minutes)
 - 4) Ethanol 95% (2 minutes)
3. Rinse sections in distilled water for 5 minutes.
4. Cover the sections with 0.5-1 ml of Silver Nitrate Solution and expose them to bright sunlight through a clear glass window for 20 minutes. Alternatively, place the sections under a 60–100-watt light bulb for 1 hour, or until the calcium turns black.

Note: If the stain is weak or fades, this indicates insufficient light exposure. In such cases, extend the incubation time as needed, up to several hours.

5. Rinse the sections thoroughly with multiple changes of distilled water.
6. To remove any unreacted silver, cover the sections with 0.5-1 ml of sodium thiosulfate and let it sit for 5 minutes. Rinse thoroughly in distilled water.
7. Counterstain with nuclear fast red for 5 minutes.
8. Rinse in distilled water.
9. Dehydrate with 2 changes of 95% Ethanol and 2 changes of 100% Ethanol (2 minutes per change).
10. Clear with 3 changes of xylene (5 minutes per change)
11. Mount coverslip onto glass slide with Permount or some other suitable organic mounting medium

Notes

1. UV light typically produces a stronger reaction, resulting in black calcium deposits. Exposure to a regular 60–100-watt light bulb often yields a brown-black coloration.
2. When confirmation is required, perform a negative control by treating a test slide with 10% formic acid for 10 minutes prior to Step 3. The treated slide should show no staining if the deposits are calcium.

Expected Results

- Calcium salts-----Black or brown black
- Nuclei-----Red
- Cytoplasm-----Pink

Positive Controls

- 16–18-day mouse embryo
- Calcium-containing tissues
- Undecalcified bone

References

1. Sheehan D, Hrapchak B, Theory and Practice of Histotechnology, 2nd Ed, 1980, pp 226-227, Battelle Press, Ohil.
2. R Marcos R, et al. Cytochemical detection of calcium in a case of calcinosis circumscripta in a dog. Vet Clin Pathol. 2006; 35(2):239-42.
3. Symonds DA, Use of the von Kossa Stain in Identifying Occult Calcifications in Breast Biopsies. American Journal of Clinical Pathology, 1990, 94 (1) 44–48,

Note: This product is intended for research purposes only. This product is not intended to be used for therapeutic or diagnostic purposes in humans or animals.

Precautions: Handle with care. Avoid contact with eyes, skin and clothing. Do not ingest. Wear gloves.